turning ideas into solutions



MB200 IGMS & MB300 IGMS INERT GAS MANAGEMENT SYSTEMS



- The market standard in high efficiency gas purifiers
- Introduced to the market in 2001 now in its 5th generation
- Developed specifically for the demands of the OLED industry
- Unmatched safety, automation and capacity characteristics
- For enclosures up to 4m³ (MB200 IGMS) and 10m³ (MB300 IGMS)
- Attainable purity of oxygen (O_2) and moisture (H_2O) < 1 ppm

www.mbraun.com

IGMS GAS PURIFICATION

MBRAUN

Working under pure atmosphere with oxygen and moisture levels below 1ppm is the prerequisite of processing high sensitive materials in the OLED industry.

The MBraun fully automatic IGMS gas purification systems are the worldwide recognized industrial standard in OLED production. A frequency controlled blower and two redundant reactors are used for continuous removal of oxygen and moisture during gas circulation.

The working pressure inside the inert gas enclosure is automatically controlled within a programmable range in over pressure as well as in under pressure. All process relevant parameters like oxygen, moisture and pressure are monitored by MBraun's proprietary designed sensors. In contrast to other commonly available gas purifiers the IGMS features a full control of all connected utilities like working gas, control gas and regeneration gas.

An automatic leak test (ALAS-test) can be performed to verify the leak tightness of the complete inert gas enclosure system.

The automatically controlled purge function and supervised change over to circulation guarantees the highest operational standard for the start up as well as during processing.

Optionally the MBraun IGMS can be equipped with a redundant solvent vapor removal system (LMF) to separate hydrocarbons from the inert gas atmosphere. The remaining capacity of the filters is detected by a special solvent sensor.

MB200 IGMS & MB300 IGMS GAS PURIFIER-E2011/04 by Daniel Kacerovsky • © M.

Further upgrades like multi-box-operation, vacuum interlock for working in combination with vacuum chambers, extended PLC control units (e.g. Mitsubishi Q-Series) or customized communication interfaces promote the MBraun IGMS to an unmatched allrounder in the inert gas technology.

MB200 IGMS & MB300 IGMS Overview	MB200 IGMS	MB200 IGMS-LMF	MB300 IGMS	MB300 IGMS-LMF
Performance				10
Max. enclosure volume [m ³]	4	4	10	10
Blower capacity [m ³ /h]	84	84	120	120
Purity O ₂ [ppm]	<1	<1	<1	<1
Purity H ₂ O [ppm]	<1	<1	<1	<1
Capacity O ₂ [NI]	100	100	275	275
Capacity H ₂ 0 [g] ⁽¹⁾	2.400	2.400	7.000	7.000
Capacity LMF [g] (2)	-	1.600	-	4.000
Sensors				
O ₂ ppm	•	•	•	•
0 ₂ %	•	•	•	•
H ₂ O ppm	•	•	•	•
Pressure sensor	•	•	•	•
Solvent sensor	-	•	-	•
Utility monitoring Features	•	•	•	•
Automatic purge [l/min.]	200	200	500	500
ALAS leak test	•	٠	٠	•
Trend diagram H ₂ O/O ₂ ppm	•	٠	٠	•
Trend diagram box pressure	•	٠	•	•
Automatic regeneration	•	•	•	•
Mitsubishi control (FX-series)	•	•	•	•
Multi color touch panel	٠	٠	•	•
Dry vacuum pump	•	•	•	•
Options				
Integrated 2-Box Operation	0	0	0	O ⁽³⁾
Integrated 3-Box Operation	0	0	0	O ⁽³⁾
Mitsubishi Q-Series control	0	0	0	0
MB-QP (Quick purge)	0	0	0	0
MB-APU (Air purge)	0	0	0	0
MB-SAPU	-	-	0	0
Vacuum interlock VCR fittings	0	0	0	0
VOR IIIIIIIgs	0	0	0	0

M.Braun Inertgas-Systeme GmbH (Headquarters)

Dieselstr. 31 · 85748 Garching · Germany Phone: +49 89 32669-0 · Fax: +49 89 32669-105 Web: www.mbraun.com · E-mail: info@mbraun.de

M.Braun Incorporated

14 Marin Way · Stratham, NH · 03885 · USA Phone: +1 603 773-9333 · Fax: +1 603 773-0008 Web: www.mbraun.com · E-mail: info@mbraunusa.com turning ideas into solutions



MB600 IGMS INERT GAS MANAGEMENT SYSTEMS



- The market standard in high efficiency gas purifiers
- Introduced to the market in 2008 now in its 2nd generation
- Developed specifically for the demands of OLED manufacturing
- Unmatched safety, automation and capacity characteristics
- For industrial enclosures up to 20m³
- Attainable purity of oxygen (O2) and moisture (H2O) < 1 ppm

www.mbraun.com

IGMS GAS PURIFICATION

Inn) **Braun**

The rise of OLED technology from R&D to large scale manufacturing defined the demands of enhanced inert gas equipment. Addressing this marked driven need the MBRAUN gas purification system MB600 IGMS was developed with the goal to purify industrial inert gas enclosures of up to 20m³ whilst ensuring the well-established MBRAUN IGMS concept.

Utilizing MBRAUN's standardized components of the MB300 IGMS such as filter units, blowers, heat exchangers and control units aligned with the proven IGMS operating software ensures a high reliability and a stable performance.

All process relevant parameters like oxygen, moisture and pressure are monitored by MBraun's proprietarily designed sensors. In contrast to other commonly available gas purifiers the IGMS features a full control of all connected utilities like working gas, control gas and regeneration gas.

Functions like the automatic leak test (ALAS) and the automatic purge function offer a maximum of process reliability.

For restricted spaces a compact version is available - the MB600 IGMS-RH. Its reduced overall height and a reinforced frame with castors the MB600 IGMS-RH allows easy move in and uncomplicated installation.

With its capacity and flexibility offered by additional options the MB600 IGMS is the system of choice to fulfill the requirements of industrial OLED production.

MB600 IGMS	SM	S
&	0 16	0 10
MB300 IGMS Overview	MB300 IGMS	MB600 IGMS
Dorformanaa	E	-
Performance Max. enclosure volume [m ³]	10	20
Blower capacity [m ³ /h]	120	200
Purity O ₂ [ppm]	<1	<1
Purity H ₂ O [ppm]	<1	<1
Capacity O ₂ [NI]	275	550
Capacity H ₂ 0 [g] ⁽¹⁾	7,000	14,000
Capacity LMF [g] (2)	-	-
Sensors		
O ₂ ppm	٠	•
O ₂ %	•	•
H ₂ O ppm	•	•
Pressure sensor	٠	•
Solvent sensor	•	-
Utility monitoring	•	•
Features Automatic purge [l/min.]	500	500
Automatic purge [//min.j	•	•
Trend diagram H_2O/O_2 ppm	•	•
Trend diagram box pressure	•	•
Automatic regeneration	•	•
Mitsubishi control (FX-series)	•	•
Multi color touch panel	•	•
Dry vacuum pump	•	•
Options		
Integrated 2-Box Operation	0	-
Integrated 3-Box Operation	0	-
Mitsubishi Q-Series control	0	0
MB-QP (Quick purge)	0	0
MB-APU (Air purge)	0	0
MB-SAPU	0	-
Vacuum interlock	0	0
VCR fittings (1) per mfg specs (2) e.g. Toulene (3) customized	0	o vailable [-] option not available

M. Braun Inertgas-Systeme GmbH (Headquarters)

Dieselstr. 31 · 85748 Garching · Germany Phone: +49 89 32669-0 · Fax: +49 89 32669-105 Web: www.mbraun.com · E-mail: info@mbraun.de

M. Braun Incorporated

14 Marin Way · Stratham, NH · 03885 · USA Phone: +1 603 773-9333 · Fax: +1 603 773-0008 Web: www.mbraun.com · E-mail: info@mbraunusa.com